



INVICTA
EDUCATION

SCIENCE
UNDERSTANDING NATURE

POOTER

PRODUCT GUIDE

IP 040359

THIS PRODUCT GUIDE MAY BE PHOTOCOPIED FOR CLASS USE ONLY

ASSEMBLY

Fit the lid onto the viewing chamber. Insert the green tube into the hole with the mesh covering and insert the clear tube into the hole without the mesh covering.


The Pooter is now ready to use.

CLEANING

We recommend that the green tube is only used by one child (you could label them) and is cleaned following each use. The tubes can be washed in soapy water or sterilised using a dilute solution of sterilising liquid (e.g. Milton). The viewing chamber, the lid and the clear tube can also be cleaned in the same way.

HEALTH & SAFETY

- Do not use the Pooter to collect specimens from water.
- Do not allow children to suck insects and small organisms from areas which are very dusty where they are likely to inhale particles into their lungs.
- Do not allow children to suck insects and small organisms from areas where there are likely to be spores as this might cause respiratory problems.
- Always wash or sterilise the Pooter after use.

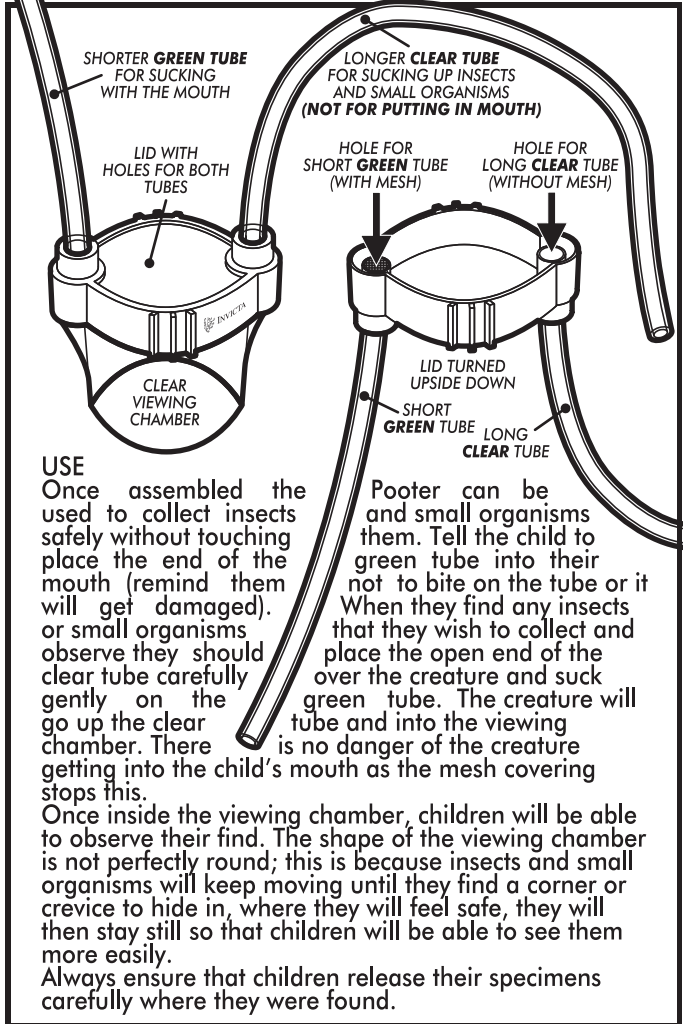

USE
Once assembled the Pooter can be used to collect insects and small organisms safely without touching them. Tell the child to place the end of the green tube into their mouth (remind them not to bite on the tube or it will get damaged). When they find any insects that they wish to collect and observe they should place the open end of the clear tube carefully over the creature and suck gently on the green tube. The creature will go up the clear tube and into the viewing chamber. There is no danger of the creature getting into the child's mouth as the mesh covering stops this.

Once inside the viewing chamber, children will be able to observe their find. The shape of the viewing chamber is not perfectly round; this is because insects and small organisms will keep moving until they find a corner or crevice to hide in, where they will feel safe, they will then stay still so that children will be able to see them more easily.

Always ensure that children release their specimens carefully where they were found.

Pooter can be used to collect insects and small organisms safely without touching them. Tell the child to place the end of the green tube into their mouth (remind them not to bite on the tube or it will get damaged). When they find any insects that they wish to collect and observe they should place the open end of the clear tube carefully over the creature and suck gently on the green tube. The creature will go up the clear tube and into the viewing chamber. There is no danger of the creature getting into the child's mouth as the mesh covering stops this.

Once inside the viewing chamber, children will be able to observe their find. The shape of the viewing chamber is not perfectly round; this is because insects and small organisms will keep moving until they find a corner or crevice to hide in, where they will feel safe, they will then stay still so that children will be able to see them more easily.

Always ensure that children release their specimens carefully where they were found.



WARNING: NOT SUITABLE FOR CHILDREN UNDER 36 MONTHS BECAUSE SMALL PARTS MAY CAUSE A CHOKING HAZARD. TO BE USED ONLY UNDER ADULT SUPERVISION. Please retain the information on this pack for future reference. We reserve the right to alter designs and specifications (including colours and materials) if and when such changes are unavoidable. This product conforms to the safety requirements of EN71, ASTM, 16 CFR, CCC and The Canadian Hazardous Products (Toys) Regulations. **MADE IN CHINA**

